



SCAN QR CODE
FOR OTHER
LANGUAGES

OPERATING INSTRUCTIONS

EN

Translation of the Original

MVP 030-3 DC

Diaphragm Pump

Dear Customer,

Thank you for choosing a Pfeiffer Vacuum product. Your new diaphragm pump is designed to support you with its performance, perfect operation and without impacting your individual application. The name Pfeiffer Vacuum stands for high-quality vacuum technology, a comprehensive and complete range of top-quality products and first-class service. From this extensive, practical experience we have gained a large volume of information that can contribute to efficient deployment and to your personal safety.

In the knowledge that our product must avoid consuming work output, we trust that our product can offer you a solution that supports you in the effective and trouble-free implementation of your individual application.

Please read these operating instructions before putting your product into operation for the first time. If you have any questions or suggestions, please feel free to contact info@pfeiffer-vacuum.de.

Further operating instructions from Pfeiffer Vacuum can be found in the [Download Center](#) on our website.

Disclaimer of liability

These operating instructions describe all models and variants of your product. Note that your product may not be equipped with all features described in this document. Pfeiffer Vacuum constantly adapts its products to the latest state of the art without prior notice. Please take into account that online operating instructions can deviate from the printed operating instructions supplied with your product.

Furthermore, Pfeiffer Vacuum assumes no responsibility or liability for damage resulting from the use of the product that contradicts its proper use or is explicitly defined as foreseeable misuse.

Copyright

This document is the intellectual property of Pfeiffer Vacuum and all contents of this document are protected by copyright. They may not be copied, altered, reproduced or published without the prior written permission of Pfeiffer Vacuum.

We reserve the right to make changes to the technical data and information in this document.

Table of contents

1	About this manual	7
1.1	Validity	7
1.1.1	Applicable documents	7
1.1.2	Variants	7
1.2	Target group	7
1.3	Conventions	7
1.3.1	Instructions in the text	7
1.3.2	Pictographs	8
1.3.3	Stickers on the product	8
1.3.4	Abbreviations	9
2	Safety	10
2.1	General safety information	10
2.2	Safety instructions	10
2.3	Safety precautions	12
2.4	Limits of use of the product	13
2.5	Proper use	13
2.6	Foreseeable improper use	13
2.7	Personnel qualification	14
2.7.1	Ensuring personnel qualification	14
2.7.2	Personnel qualification for maintenance and repair	15
2.7.3	Advanced training with Pfeiffer Vacuum	15
3	Product description	16
3.1	Function	16
3.1.1	Actuator	16
3.1.2	Pumping system	16
3.1.3	Cooling	16
3.2	Connections	16
3.3	Identifying the product	17
3.4	Scope of delivery	17
4	Transportation and Storage	18
4.1	Transporting vacuum pump	18
4.2	Storing vacuum pump	18
5	Installation	19
5.1	Installing the vacuum pump	19
5.2	Connecting vacuum side	19
5.3	Connect exhaust side	20
5.4	Establishing electric connection	20
6	Interfaces	23
6.1	"Remote" interface	23
6.1.1	Voltage supply	24
6.1.2	Inputs	24
6.1.3	Outputs	24
6.1.4	RS-485	25
6.2	Interface RS-485	25
6.3	Cross-linking via RS-485 connection	25
6.4	Connection options via interface RS-485	26
6.5	Pfeiffer Vacuum protocol for RS-485 interface	27
6.5.1	Telegram frame	27
6.5.2	Telegram description	27
6.5.3	Telegram example 1	27
6.5.4	Telegram example 2	28
6.5.5	Data types	28

7	Parameter set	29
7.1	General	29
7.2	Control commands	29
7.3	Status requests	30
7.4	Reference value inputs	31
8	Operation	32
8.1	Commissioning vacuum pump	32
8.2	Switching on the vacuum pump	32
8.3	Configuring connections with Pfeiffer Vacuum parameter set	33
8.3.1	Configuring the digital outputs	33
8.3.2	Selecting the interfaces	33
8.4	Operating modes	33
8.4.1	Rotation speed setting operation	34
8.4.2	Standby mode	34
8.4.3	Normal operation	34
8.5	Pumping condensable vapors	35
8.6	Operation monitoring	36
8.6.1	Operating mode display via LED	36
8.6.2	Temperature monitoring	36
8.7	Switching off the vacuum pump	36
9	Maintenance	37
9.1	Maintenance instructions	37
9.2	Checklist for inspection and maintenance	38
9.3	Dismantling / installing the handle	38
9.4	Change the diaphragms and valves	39
9.4.1	Dismantle the pump head and valves	39
9.4.2	Dismantling the diaphragms	40
9.4.3	Installing the diaphragm	41
9.4.4	Installing the pump head and valves	42
9.5	Checking final pressure	42
10	Decommissioning	44
11	Recycling and disposal	45
11.1	General disposal information	45
11.2	Dispose of diaphragm pumps	45
12	Malfunctions	46
12.1	General	46
12.2	Error codes	47
13	Service solutions by Pfeiffer Vacuum	48
14	Spare part packages	50
15	Accessories	52
15.1	Accessory information	52
15.2	Ordering accessories	52
16	Technical data and dimensions	53
16.1	General	53
16.2	Technical data	53
16.3	Substances in contact with media	54
16.4	Dimensions	55
	EC Declaration of Conformity	56
	UK Declaration of Conformity	57

List of tables

Tbl. 1:	Stickers on the product	8
Tbl. 2:	Abbreviations used in this document	9
Tbl. 3:	Limits of use of the product	13
Tbl. 4:	Intake temperature limits of use	13
Tbl. 5:	Connection description of the electronic drive unit	16
Tbl. 6:	Pin assignment of the D-Sub socket, 15-pin	23
Tbl. 7:	Output DO1/pin 8	24
Tbl. 8:	Output DO2/pin 9	24
Tbl. 9:	Features of the RS-485 interface	25
Tbl. 10:	Explanation and meaning of the parameters	29
Tbl. 11:	Parameter set Control commands	30
Tbl. 12:	Parameter set Status requests	30
Tbl. 13:	Parameter set Reference value inputs	31
Tbl. 14:	Factory setting of the electronic drive unit for diaphragm pumps when delivered	32
Tbl. 15:	Configure parameters [P:019] and [P:024]	33
Tbl. 16:	Parameters [P:060] and [P:061]	33
Tbl. 17:	Behavior and meaning of the LED of the electronic drive unit	36
Tbl. 18:	Typical service life with normal use	37
Tbl. 19:	Maintenance intervals	38
Tbl. 20:	Troubleshooting on diaphragm pumps	47
Tbl. 21:	Error messages for vacuum pump	47
Tbl. 22:	Warning messages for vacuum pump	47
Tbl. 23:	Spare part packages	50
Tbl. 24:	Spart part package flushing gas nozzle	51
Tbl. 25:	Accessories	52
Tbl. 26:	Conversion table: Pressure units	53
Tbl. 27:	Conversion table: Units for gas throughput	53
Tbl. 28:	Technical data, MVP 030-3 DC	54
Tbl. 29:	Materials that make contact with the process media	54

List of figures

Fig. 1:	Position of the stickers on the product	8
Fig. 2:	Diaphragm pump design	16
Fig. 3:	Minimum distances	19
Fig. 4:	Connection with turbopump, power supply pack and controller	21
Fig. 5:	Connection to control unit with integrated power supply pack	21
Fig. 6:	Pins of the D-Sub socket, 15-pin	23
Fig. 7:	Cross-linking via RS-485 interface	26
Fig. 8:	Connection options via interface RS-485	26
Fig. 9:	Flushing gas nozzle with filter	35
Fig. 10:	Handle	39
Fig. 11:	Pump head and valves	40
Fig. 12:	Replacing the diaphragms	41
Fig. 13:	Correct mounting orientation of the valves in the head cover	42
Fig. 14:	Spare parts	50
Fig. 15:	Dimensions MVP 030-3 DC	55

1 About this manual



IMPORTANT

Read carefully before use.
Keep the manual for future consultation.

1.1 Validity

These operating instructions are a customer document of Pfeiffer Vacuum. The operating instructions describe the functions of the named product and provide the most important information for the safe use of the device. The description is written in accordance with the valid directives. The information in these operating instructions refers to the product's current development status. The document shall remain valid provided that the customer does not make any changes to the product.

1.1.1 Applicable documents

Designation	Document
Declaration of conformity	A component of these operating instructions

1.1.2 Variants

These instructions apply to diaphragm pumps of the DC series:

- MVP 030-3 DC as standard version
- MVP 030-3 DC in the version for the integration in a pumping station (without handle and rubber feet)

1.2 Target group

These operating instructions are aimed at all persons performing the following activities on the product:

- Transportation
- Setup (Installation)
- Usage and operation
- Decommissioning
- Maintenance and cleaning
- Storage or disposal

The work described in this document is only permitted to be performed by persons with the appropriate technical qualifications (expert personnel) or who have received the relevant training from Pfeiffer Vacuum.

1.3 Conventions

1.3.1 Instructions in the text

Usage instructions in the document follow a general structure that is complete in itself. The required action is indicated by an individual step or multi-part action steps.

Individual action step

A horizontal, solid triangle indicates the only step in an action.

- This is an individual action step.

Sequence of multi-part action steps

The numerical list indicates an action with multiple necessary steps.

1. Step 1
2. Step 2
3. ...

1.3.2 Pictographs

Pictographs used in the document indicate useful information.



Note



Tip

1.3.3 Stickers on the product

This section describes all the stickers on the product along with their meaning.

	Rating plate Rating plate of the diaphragm pump
	Warning of hot surfaces The sticker warns of hot surfaces during operation or immediately after switching off the vacuum pump.
	General warning signs The sticker provides a general warning of possible dangers.
	Closure seal The diaphragm heads are sealed at the factory. Damaging or removing a closure seal results in loss of the warranty.

Tbl. 1: Stickers on the product

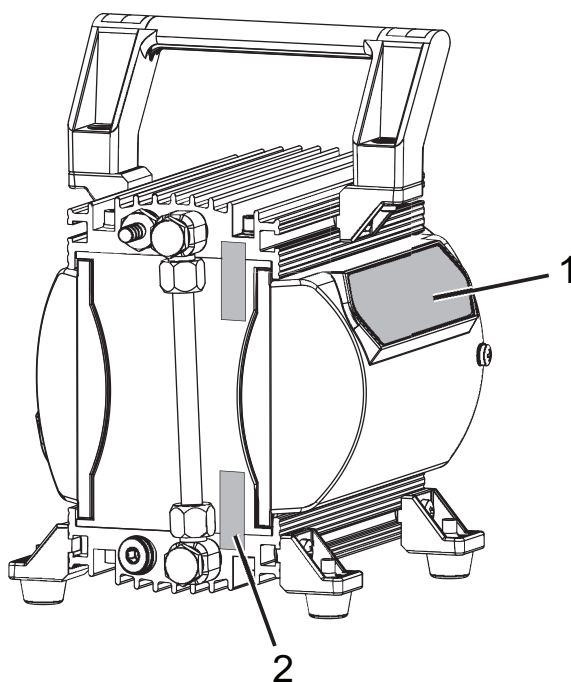


Fig. 1: Position of the stickers on the product

1 Rating plate of the diaphragm pump
Incl. the 2 warning signs

2 Warranty seal (2×)

1.3.4 Abbreviations

Abbreviation	Meaning in this document
DC	Direct current
DN	Nominal diameter (diamètre nominal)
EMC	Electromagnetic compatibility
f	Rotation speed value of a vacuum pump (frequency, in rpm or Hz)
FKM	Fluorinated rubber
GND	Ground
LED	Light emitting diode
MVP	Diaphragm vacuum pump
MSL	Mean sea level
[P:xxx]	Electronic drive unit control parameters. Printed in bold as a three-digit number in square brackets. Frequently displayed in conjunction with a short description. Example: [P:312] software version
PE	Protective earth (earthed conductor)
PWM	Pulse width modulation (Type of modulation where one technical variable (e.g. electrical voltage) changes between two values.)
T	Temperature (in °C), property abbreviation of the vacuum pump
TC	Turbopump electronic drive unit (turbo controller)
TPS	Voltage supply (turbo power supply)

Tbl. 2: Abbreviations used in this document

2 Safety

2.1 General safety information

The following 4 risk levels and 1 information level are taken into account in this document.

DANGER

Immediately pending danger

Indicates an immediately pending danger that will result in death or serious injury if not observed.

- Instructions to avoid the danger situation

WARNING

Potential pending danger

Indicates a pending danger that could result in death or serious injury if not observed.

- Instructions to avoid the danger situation

CAUTION

Potential pending danger

Indicates a pending danger that could result in minor injuries if not observed.

- Instructions to avoid the danger situation

NOTICE

Danger of damage to property

Is used to highlight actions that are not associated with personal injury.

- Instructions to avoid damage to property



Notes, tips or examples indicate important information about the product or about this document.

2.2 Safety instructions

All safety instructions in this document are based on the results of the risk assessment carried out in accordance with Machinery Directive 2006/42/EC Annex I and EN ISO 12100 Section 5. As far as applicable, all unit life cycle phases have been considered.

Risks during transport

WARNING

Danger of serious injury due to falling objects

Due to falling objects there is a risk of injuries to limbs through to broken bones.

- Take particular care and pay special attention when transporting products manually.
- Do not stack the products.
- Wear protective equipment, e.g. safety shoes.

Risks during installation

DANGER

Danger to life from electric shock

Applying voltages that exceed the specified safety extra-low voltage (according to IEC 60449 and VDE 0100) will result in the destruction of the insulation. There is a danger to life from electric shock at the communication interfaces.

- Connect only suitable devices to the bus system.

⚠ CAUTION**Danger of injury from bursting as a result of high pressure in the exhaust duct**

Faulty or inadequate exhaust ducts lead to dangerous situations, e.g., increased exhaust pressure. There is a danger of bursting. Injuries caused by flying fragments, the escaping of high pressure, and damage to the unit cannot be excluded.

- ▶ Route the exhaust duct without shut-off units.
- ▶ Observe the permissible pressures and pressure differentials for the product.
- ▶ Check the proper function of the exhaust duct on a regular basis.

Risks during operation**⚠ WARNING****Danger of poisoning due to toxic process media escaping from the exhaust duct.**

During operation with no exhaust duct, the vacuum pump allows exhaust gases and vapors to escape freely into the air. There is a risk of injury and fatalities due to poisoning in processes with toxic process media.

- ▶ Observe the pertinent regulations for handling toxic process media.
- ▶ Safely purge toxic process media via an exhaust duct.
- ▶ Use appropriate filter equipment to separate toxic process media.

⚠ WARNING**Risk of explosion from reactive, explosive or other hazardous gas / air mixtures.**

Uncontrolled inlet of gas at the flushing gas nozzle leads to sparks generated mechanically following diaphragm rupture, hot surfaces or to explosive gas / air mixtures in the vacuum system caused by static electricity.

- ▶ If necessary, use inert gas for supplying the flushing gas in order to avoid a potential ignition.

⚠ CAUTION**Danger of burns on hot surfaces**

Depending on the operating and ambient conditions, the surface temperature of the vacuum pump can increase to above 70 °C.

- ▶ Provide suitable touch protection.

⚠ CAUTION**Risk of poisoning due to toxic process gases escaping**

For intake pressures > 500 hPa, process gas may escape from the flushing gas nozzle.

- ▶ If necessary, use an extraction line.

Risks during maintenance, decommissioning and in the event of malfunctions**⚠ WARNING****Health hazard through poisoning from toxic contaminated components or devices**

Toxic process media result in contamination of devices or parts of them. During maintenance work, there is a risk to health from contact with these poisonous substances. Illegal disposal of toxic substances causes environmental damage.

- ▶ Take suitable safety precautions and prevent health hazards or environmental pollution by toxic process media.
- ▶ Decontaminate affected parts before carrying out maintenance work.
- ▶ Wear protective equipment.

CAUTION

Danger of injury from moving parts

After a power failure or a standstill as a result of overheating, the motor restarts automatically. There is a risk of injury to fingers and hands if they enter the operating range of rotating parts.

- ▶ Safely disconnect motor from the mains.
- ▶ Secure the motor against reactivation.
- ▶ Dismantle the vacuum pump for inspection, away from the system if necessary.

CAUTION

Danger of burns on hot surfaces

In the event of a fault, the surface temperature of the vacuum pump can increase to above 105°C.

- ▶ Allow the vacuum pump to cool down before carrying out any work.
- ▶ Wear personal protective equipment if necessary.

Risks during disposal

WARNING

Health hazard through poisoning from toxic contaminated components or devices

Toxic process media result in contamination of devices or parts of them. During maintenance work, there is a risk to health from contact with these poisonous substances. Illegal disposal of toxic substances causes environmental damage.

- ▶ Take suitable safety precautions and prevent health hazards or environmental pollution by toxic process media.
- ▶ Decontaminate affected parts before carrying out maintenance work.
- ▶ Wear protective equipment.

2.3 Safety precautions

DANGER

Danger to life from electric shock

Power supply packs that are not specified or are not approved will lead to severe injury to death.

- ▶ Make sure that the power supply pack meets the requirements for double isolation between mains input voltage and output voltage, in accordance with IEC 61010-1 IEC 60950-1 and IEC 62368-1.
- ▶ Make sure that the power supply pack meets the requirements in accordance with IEC 61010-1 IEC 60950-1 and IEC 62368-1.
- ▶ Where possible, use original power supply packs or only power supply packs that correspond with the applicable safety regulations.



Duty to provide information on potential dangers

The product holder or user is obliged to make all operating personnel aware of dangers posed by this product.

Every person who is involved in the installation, operation or maintenance of the product must read, understand and adhere to the safety-related parts of this document.



Infringement of conformity due to modifications to the product

The Declaration of Conformity from the manufacturer is no longer valid if the operator changes the original product or installs additional equipment.

- Following the installation into a system, the operator is required to check and re-evaluate the conformity of the overall system in the context of the relevant European Directives, before commissioning that system.

General safety precautions

- ▶ Prevent process gases and their reaction and ancillary products that are harmful to health and the environmental from escaping, or dispose of these according to the applicable regulations.
- ▶ Avoid contact with process gases that are harmful to health and if required, wear personal protective clothing.
- ▶ Check the media compatibility of parts carrying media.
- ▶ Observe all applicable safety and accident prevention regulations.
- ▶ Check that all safety measures are observed at regular intervals.
- ▶ Do not expose body parts to the vacuum.
- ▶ Never disconnect plug connections during operation.
- ▶ Never fill or operate the unit with cleaning agents or cleaning agent residues.
- ▶ Do not carry out your own conversions or modifications on the unit.
- ▶ Observe the unit protection degree prior to installation or operation in other environments.
- ▶ Provide suitable touch protection, if the surface temperature exceeds 70°C.
- ▶ Keep lines and cables away from hot surfaces.
- ▶ Before returning the vacuum pump, observe the notes in the chapter Service solutions.

2.4 Limits of use of the product

Installation location	Interiors, protected against the accumulation of dust and weather influences, in a non-explosive, dry environment
Protection degree IEC 60529	IP 20
Protection degree UL 50E	NEMA type 1
Installation altitude	max. 2000 m above sea level ¹⁾
Ambient temperature	+10 °C to +40 °C
Relative air humidity	80% at T ≤ 31 °C, to max. 50% at T ≤ 40 °C
Degree of contamination	2

Tbl. 3: Limits of use of the product

Operations condition, status	Intake pressure	Permissible intake temperature of medium to be pumped
Continuous operation	> 100 hPa (high gas load)	+10 °C to +40 °C
Continuous operation	< 100 hPa (low gas load)	0 °C to +60 °C
short-term (< 5 minutes)	< 100 hPa (low gas load)	-10 °C to +80 °C

Tbl. 4: Intake temperature limits of use

2.5 Proper use

- ▶ Use the vacuum pump for vacuum generation only.
- ▶ Adhere to the installation, commissioning, operating, and maintenance instructions.
- ▶ Use only accessory parts recommended by Pfeiffer Vacuum.

2.6 Foreseeable improper use

Improper use of the product invalidates all warranty and liability claims. Any use that is counter to the purpose of the product, whether intentional or unintentional, is regarded as improper use; in particular:

- Pumping of corrosive media
- Pumping explosive media
- Pumping radioactive or volatile media
- Pumping of gases that contain contamination such as particles, dust, or condensate
- Pumping of fluids

¹⁾ at an installation location higher than 1000 meters above sea level, there is a risk of insufficient cooling. If necessary, take measures according to Directive DIN EN 61010.

- Pumping FKM-solubilizing media
- Pumping vapors from combustible liquids
- Pumping pressurized media (> atmospheric pressure)
- Pumping media that can condense or cause adhering deposits to form in the suction chamber
- Using the vacuum pump outside the specified range of application
- Using the vacuum pump below ground
- Using the vacuum pump to generate pressure
- Use of the vacuum pump in systems in which sporadic loads and vibrations or periodic forces act on the unit
- Using the vacuum pump in strong electrical, magnetic, or electromagnetic fields
- Connecting to vacuum pumps and units that are not designed for this purpose according to their operating instructions
- Connecting to units with exposed live parts
- Connecting to sockets without earthing contact
- Use of pipes to lift the vacuum pump
- Use of accessories or spare parts not listed in these instructions
- Using the vacuum pump as a climbing aid
- Using the connection lines between the diaphragm heads as carrying handles
- Using the pump in ambient conditions that do not meet the specified IP protection degree limits

2.7 Personnel qualification

The work described in this document may only be carried out by persons who have appropriate professional qualifications and the necessary experience or who have completed the necessary training as provided by Pfeiffer Vacuum.

Training people

1. Train the technical personnel on the product.
2. Only let personnel to be trained work with and on the product when under the supervision of trained personnel.
3. Only allow trained technical personnel to work with the product.
4. Before starting work, make sure that the commissioned personnel have read and understood these operating instructions and all applicable documents, in particular the safety, maintenance and repair information.

2.7.1 Ensuring personnel qualification

Specialist for mechanical work

Only a trained specialist may carry out mechanical work. Within the meaning of this document, specialists are people responsible for construction, mechanical installation, troubleshooting and maintenance of the product, and who have the following qualifications:

- Qualification in the mechanical field in accordance with nationally applicable regulations
- Knowledge of this documentation

Specialist for electrotechnical work

Only a trained electrician may carry out electrical engineering work. Within the meaning of this document, electricians are people responsible for electrical installation, commissioning, troubleshooting, and maintenance of the product, and who have the following qualifications:

- Qualification in the electrical engineering field in accordance with nationally applicable regulations
- Knowledge of this documentation

In addition, these individuals must be familiar with applicable safety regulations and laws, as well as the other standards, guidelines, and laws referred to in this documentation. The above individuals must have an explicitly granted operational authorization to commission, program, configure, mark, and earth devices, systems, and circuits in accordance with safety technology standards.

Trained individuals

Only adequately trained individuals may carry out all works in other transport, storage, operation and disposal fields. Such training must ensure that individuals are capable of carrying out the required activities and work steps safely and properly.

2.7.2 Personnel qualification for maintenance and repair



Advanced training courses

Pfeiffer Vacuum offers advanced training courses to maintenance levels 2 and 3.

Adequately trained individuals are:

- **Maintenance level 1**
 - Customer (trained specialist)
- **Maintenance level 2**
 - Customer with technical education
 - Pfeiffer Vacuum service technician
- **Maintenance level 3**
 - Customer with Pfeiffer Vacuum service training
 - Pfeiffer Vacuum service technician

2.7.3 Advanced training with Pfeiffer Vacuum

For optimal and trouble-free use of this product, Pfeiffer Vacuum offers a comprehensive range of courses and technical trainings.

For more information, please contact [Pfeiffer Vacuum technical training](#).

3 Product description

3.1 Function

The diaphragm pump of the MVP 030-3 DC series is a dry compressing vacuum pump with 3 pumping stages. The vacuum pump is a positive displacement pump with a periodic change in suction chamber size, produced by the movement of the diaphragm. The gas flow causes the valves to open and close automatically. The pump units are directly connected to the drive motor.

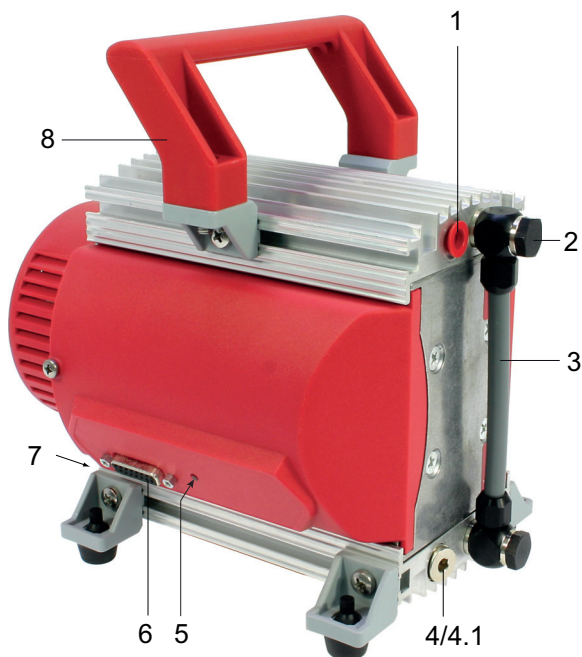


Fig. 2: Diaphragm pump design

- | | |
|------------------------------------|-------------------------|
| 1 Vacuum connection, thread 1/8" | 5 LED |
| 2 Banjo bolt | 6 Connecting socket |
| 3 Interhead connection | 7 Exhaust with silencer |
| 4 Locking screw | 8 Handle |
| 4.1 Flushing gas nozzle (optional) | |

3.1.1 Actuator

The diaphragm pump forms a compact unit with the electronic drive unit. The integrated electronic drive unit permits the control and monitoring of pumping functions via the "remote" interface using a Pfeiffer Vacuum control unit.

3.1.2 Pumping system

The diaphragm pump has 4 diaphragm heads and 3 pumping stages.

3.1.3 Cooling

The diaphragm pump is convection-cooled. The drive motor has its own fan.

3.2 Connections

Connection	Description
	remote D-sub socket with 15 pins for connection and configuration of a remote control.

Tbl. 5: Connection description of the electronic drive unit

3.3 Identifying the product

- ▶ To ensure clear identification of the product when communicating with Pfeiffer Vacuum, always keep all of the information on the rating plate to hand.
- ▶ Learn about certifications through test seals on the product or at www.certipedia.com with company ID no. 000021320.

3.4 Scope of delivery

- Vacuum pump with integrated electronic drive unit
- Silencer
- Operating instructions
- Blind plug on the vacuum connection

4 Transportation and Storage

4.1 Transporting vacuum pump

WARNING

Danger of serious injury due to falling objects

Due to falling objects there is a risk of injuries to limbs through to broken bones.

- ▶ Take particular care and pay special attention when transporting products manually.
- ▶ Do not stack the products.
- ▶ Wear protective equipment, e.g. safety shoes.



Packing

We recommend keeping the transport packaging and original protective cover.

Safe transport of the product

1. Observe the weight specified on the packaging.
2. Use a transport aid if necessary (trolley, lift truck).
3. Transport the product in its original packaging.
4. Always place the product on an adequately sized, level surface.

Transporting the vacuum pump without its packaging

1. Unpack the vacuum pump.
2. To protect the inside of the pump, leave both protective caps on the connections during transport.
3. For lifting, use the handle provided for this purpose, located on the top of the pump.
4. Lift the vacuum pump out of the transport packaging.
5. Make sure that no forces are acting on the piping system.
6. Always place the vacuum pump on an adequately sized, level surface.

4.2 Storing vacuum pump



Packing

We recommend storing the product in its original packaging.

Safe storing vacuum pump

- ▶ Seal the vacuum connection with the blind plug.
- ▶ Store the vacuum pump only in dry, dust-free rooms, within the specified ambient conditions.
- ▶ In rooms with humid or aggressive atmospheres: Hermetically seal the vacuum pump together with a drying agent in a plastic bag.

5 Installation

5.1 Installing the vacuum pump

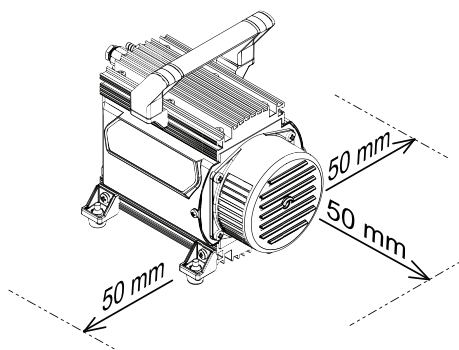


Fig. 3: Minimum distances

Procedure

1. Place the vacuum pump on a flat, horizontal surface.
2. For stationary installation, screw the vacuum pump directly to the base if necessary.
 - For this purpose, dismantle the rubber feet.
3. When installing the pump in a closed housing, ensure adequate air circulation.
 - The vacuum pump does not have a fan.
4. Make sure that minimum distances to adjacent components, e.g., walls, housings, etc., are observed.
5. Keep the specifications on the motor rating plate visible and freely accessible.

5.2 Connecting vacuum side

NOTICE

Property damage from contaminated gases

Pumping gases that contain contamination damages the vacuum pump.

- Use suitable filters or separators from the Pfeiffer Vacuum range of accessories, to protect the vacuum pump.



Installation and operation of accessories

Pfeiffer Vacuum offers a series of special, compatible accessories for its diaphragm pumps.

- Information and ordering options for approved [accessories](#) can be found online.
- Described accessories are not included in the shipment.



Preventing throttling losses

Using short vacuum lines with a large nominal diameter prevents throttling losses.



Condensate separator

Pfeiffer Vacuum recommends the installation of a condensate separator in case vapors are formed from moisture during evacuation.

Procedure

1. Remove the blind plug from the vacuum connection.
2. Establish the shortest possible connection between vacuum pump and vacuum chamber.
3. Choose a minimum vacuum line diameter equal to the nominal diameter of the vacuum connection.
4. Depending on the pump type, use PVC or metallic hoses with flange connections from the [Pfeiffer Vacuum component shop](#).
5. Connect the vacuum pump to the vacuum system using the vacuum connection.

5.3 Connect exhaust side

WARNING

Danger to life from poisoning where toxic process gases leak with no exhaust line

During normal operation, the vacuum pump expels exhaust gases and vapors freely into the air. In processes involving toxic media, there is a risk of injury and danger to life due to poisoning.

- ▶ Observe the relevant regulations for handling toxic substances.
- ▶ Safely purge toxic process gases via an exhaust line.

CAUTION

Danger of injury from bursting as a result of high pressure in the exhaust duct

Faulty or inadequate exhaust ducts lead to dangerous situations, e.g., increased exhaust pressure. There is a danger of bursting. Injuries caused by flying fragments, the escaping of high pressure, and damage to the unit cannot be excluded.

- ▶ Route the exhaust duct without shut-off units.
- ▶ Observe the permissible pressures and pressure differentials for the product.
- ▶ Check the proper function of the exhaust duct on a regular basis.

CAUTION

Health hazard from increased noise emission

Operation without silencer leads to higher noise emissions. Remaining in the close proximity of the vacuum pump for a sustained period of time may cause hearing damage.

- ▶ Install a suitable exhaust line.
- ▶ Wear hearing protection.



Condensate separator

Pfeiffer Vacuum recommends installing a condensate separator, with condensate drain at the lowest point of the exhaust line.

Procedure

1. Check the installed silencer for free passage.
2. Choose a minimum exhaust line diameter equal to the nominal diameter of the connection flange.
3. Depending on the pump type, use PVC or metallic hoses with flange connections from the [Pfeiffer Vacuum component shop](#).
4. Option: In case of higher gas throughputs, mount an exhaust line.
5. Route the piping downwards from the vacuum pump, to prevent condensate return.
6. Support or suspend the piping to the vacuum pump so that no piping system forces act on the vacuum pump.

5.4 Establishing electric connection

DANGER

Danger to life from electric shock

Power supply packs that are not specified or are not approved will lead to severe injury to death.

- ▶ Make sure that the power supply pack meets the requirements for double isolation between mains input voltage and output voltage, in accordance with IEC 61010-1 IEC 60950-1 and IEC 62368-1.
- ▶ Make sure that the power supply pack meets the requirements in accordance with IEC 61010-1 IEC 60950-1 and IEC 62368-1.
- ▶ Where possible, use original power supply packs or only power supply packs that correspond with the applicable safety regulations.

⚠ WARNING

Risk of injury due to incorrect installation

Dangerous situations may arise from unsafe or incorrect installation.

- ▶ Do not carry out your own conversions or modifications on the unit.
- ▶ Ensure the integration into an Emergency Off safety circuit.

Original power supply packs (e.g. TPS 110) and connection cables are available for the electronic drive unit supply voltage.

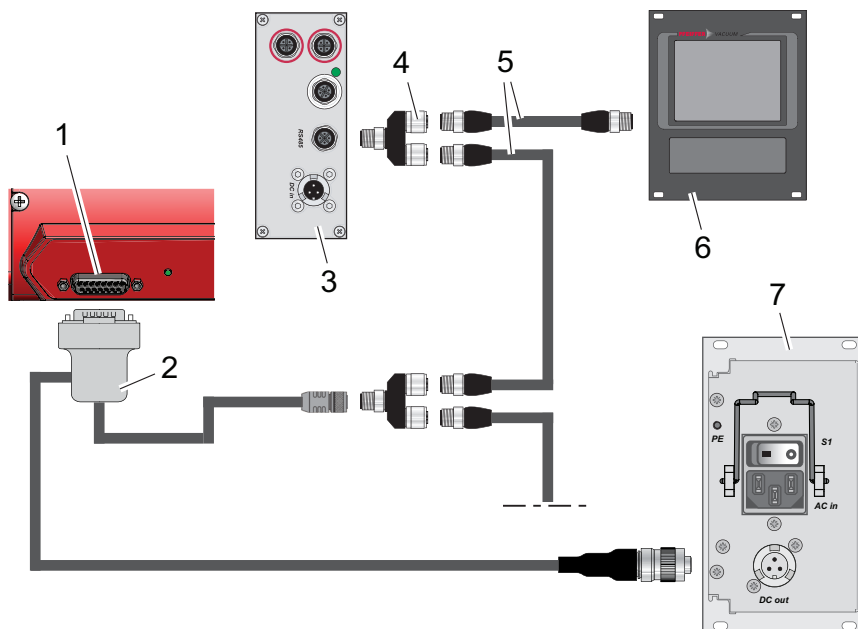


Fig. 4: Connection with turbopump, power supply pack and controller

- | | |
|---|----------------------------|
| 1 "Remote" connection on MVP electronic drive unit | 5 Connection cable M12 |
| 2 Connection cable for TPS power supply pack with RS-485 interface (order number PM 061 350 -T) | 6 OmniControl control unit |
| or
Connection cable for TC and TPS (order number PE 100 013 -T) | |
| 3 Turbopump electronic drive unit (TC) | 7 Power supply pack TPS |
| 4 Y-distributor M12 for RS-485 interface (order number P 4723 010) | |

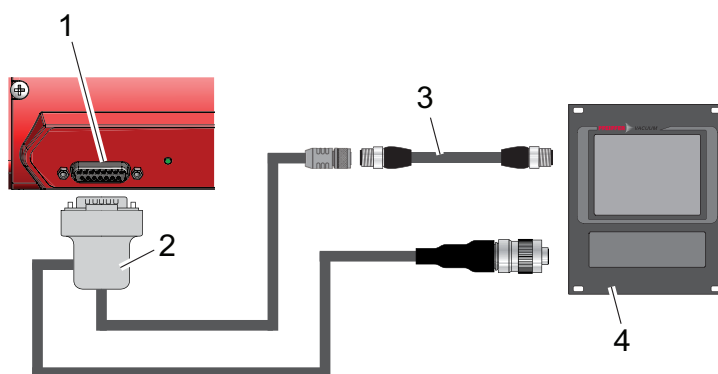


Fig. 5: Connection to control unit with integrated power supply pack

- | | |
|---|----------------------------|
| 1 "Remote" connection on MVP electronic drive unit | 3 Connection cable M12 |
| 2 Connection cable for OmniControl with RS-485 interface (order number PM 061 350 -T) | 4 OmniControl control unit |

Connecting MVP electronic drive unit

- ▶ Ensure the correct supply voltage.
- ▶ Make sure that the power supply pack main switch is off prior to connection.

- ▶ Use a suitable connection cable from the Pfeiffer Vacuum accessories range.
- ▶ Plug the 15-pin connection cable plug into the "remote" connection on the MVP electronic drive unit and secure it.
- ▶ Insert the connection cable into the connection "DC out" on the power supply pack and close the bayonet lock.
- ▶ **If you are using a Pfeiffer Vacuum control unit:** Connect the "RS-485" connector to the control unit using a suitable extension cable.

6 Interfaces

6.1 "Remote" interface

NOTICE

Property damage on the electronics

Separating all plug-and-socket connections within the bus system with voltage supply switched on may lead to the destruction of electronic components.

- ▶ Always disconnect the voltage supply before removing the connecting plug.
- ▶ After switching off the power supply pack, wait until the residual load has dispersed completely before disconnecting the plug-and-socket connection.

NOTICE

Impairment of electrical connections caused by external disruptive influences

External interference caused by electromagnetic emissions causes malfunctions.

- ▶ A screened cable should be used due to the electromagnetic compatibility (EMC).
- ▶ Connect the screening to the connector housings on both sides.

The 15-pin sub-D connection with the "remote" designation offers the possibility to operate the electronic drive unit via remote control. The following specifications are the factory settings for the electronic drive unit.

Configuring "remote" interface

- ▶ Utilize the screened plug and cable.
- ▶ Configure the inputs and outputs via the Pfeiffer Vacuum parameter set.

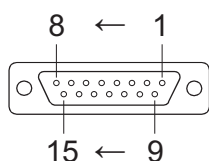


Fig. 6: Pins of the D-Sub socket, 15-pin

Pin	Function	Description, factory setting
1	+24 VDC input	Voltage supply for drive and interface
2	DI access request	V+: Control via DIs, --> GND/open: Control unlocked
3	DI1	V+: Rotation speed setting mode, --> GND/open: no rotation speed setting mode
4	DI2	Only valid for vacuum pumps with a solenoid valve; the following then applies: V+: Valve released, --> GND/open: Valve closed
5	DI pumping station	V+: Vacuum pump on, --> GND/open: Vacuum pump off
6	DI standby	V+: standby, --> GND/open: no standby
7	V+	24 V output
8	DO1	GND: Error, V+: no error
9	DO2	GND: Vacuum pump off, V+: Vacuum pump on
10, 11, 12	n.c.	-
13	RS-485	D+
14	RS-485	D-
15	Ground (GND)	Ground connection of the voltage supply; reference ground for all digital inputs and outputs

Tbl. 6: Pin assignment of the D-Sub socket, 15-pin

6.1.1 Voltage supply

Input/pin 1

The electrical connection is made using a connection cable from the Pfeiffer Vacuum accessory range or, by the customer, at pin 1 and pin 15.

+24 V DC* output/pin 7

A connection with +24 V DC to pin 7 (active high) activates inputs 2 to 6. Alternatively, they can be activated via an external PLC. "PLC High level" activates and "PLC Low level" deactivates the functions.

- PLC High level: +13 V to +33 V
- PLC Low level: -33 V to +7 V
- Ri: 7 kΩ
- $I_{\max} < 200 \text{ mA}$ (with RS-485, where present)

6.1.2 Inputs

The digital inputs switch various electronic drive unit functions. Inputs are assigned with functions ex-factory. You can configure them via the RS-485 interface and the Pfeiffer Vacuum parameter set.

Input DI remote priority/pin 2

V+: The "remote" connection has operating priority over all other control sources.

open: Remote priority "inactive"

Input DI1 rotation speed setting mode/pin 3

V+: Rotation speed setting mode "active"

open: Rotation speed setting mode "inactive"

Input DI pumping station/pin 5

V+:	Vacuum pump on Control of all components connected and malfunction acknowledgement.
open:	Vacuum pump off

Input DI stand-by/pin 6

V+:	Stand-by activated The stand-by speed can be selected in the range 30 to 100% of the nominal speed.
GND/open:	Stand-by off

6.1.3 Outputs

The digital outputs have a maximum load limit of 24 V/50 mA per output.

V+:	No error After applying the voltage supply, the digital output DO1 permanently outputs V+ meaning which means "no error".
GND	Error "Active low" signifies "Error" (common error message).

Tbl. 7: Output DO1/pin 8

V+:	Vacuum pump on "Active high" means "Vacuum pump on" and rotates at set rotation speed. Example: use the signal for the message vacuum pump "Ready for operation".
GND	Vacuum pump off

Tbl. 8: Output DO2/pin 9

6.1.4 RS-485

Connecting RS-485 via D-Sub

- Connect either a Pfeiffer Vacuum control unit or an external PC via pin 13 and pin 14 on the D-sub connector of the electronic drive unit.

6.2 Interface RS-485

DANGER

Danger to life from electric shock

Applying voltages that exceed the specified safety extra-low voltage (according to IEC 60449 and VDE 0100) will result in the destruction of the insulation. There is a danger to life from electric shock at the communication interfaces.

- Connect only suitable devices to the bus system.

The interface designated "RS-485" is intended for connecting a Pfeiffer Vacuum control unit or an external PC. The connections are galvanically safe and are isolated from the maximum supply voltage for the electronic drive unit.

Designation	Value
Serial interface	RS-485
Baud rate	9600 Baud
Data word length	8 bit
Parity	none (no parity)
Start bits	1
Stop bits	1

Tbl. 9: Features of the RS-485 interface

Connecting control units for Pfeiffer Vacuum or a PC

- Use the supplied control unit connecting cable or a cable from [Pfeiffer Vacuum's accessory range](#).
- Connect the control unit to the RS-485 interface.
- Connect a PC via the USB/RS-485 converter.

Networking as an RS-485 bus

The group address for the electronic drive unit is **902**.

1. Install the devices according to the RS-485 interface specification.
2. Make sure that all devices connected to the bus have different RS-485 device addresses **[P:797]**.
3. Connect all devices with RS-485 D+ and RS-485 D- to the bus.

6.3 Cross-linking via RS-485 connection

DANGER

Danger to life from electric shock

Applying voltages that exceed the specified safety extra-low voltage (according to IEC 60449 and VDE 0100) will result in the destruction of the insulation. There is a danger to life from electric shock at the communication interfaces.

- Connect only suitable devices to the bus system.

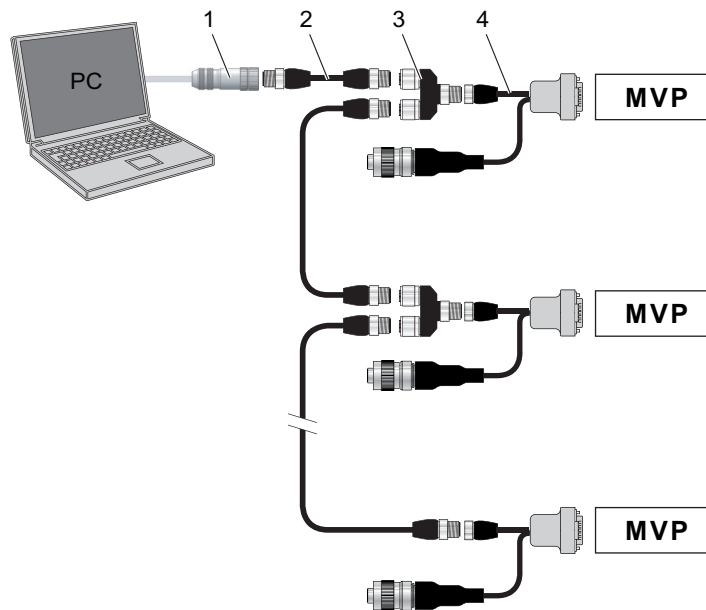


Fig. 7: Cross-linking via RS-485 interface

- | | |
|------------------------------|--------------------------------|
| 1 USB RS-485 converter | 3 Y-connector for RS-485 |
| 2 M12 to M12 extension cable | 4 Connection cable with RS-485 |

6.4 Connection options via interface RS-485

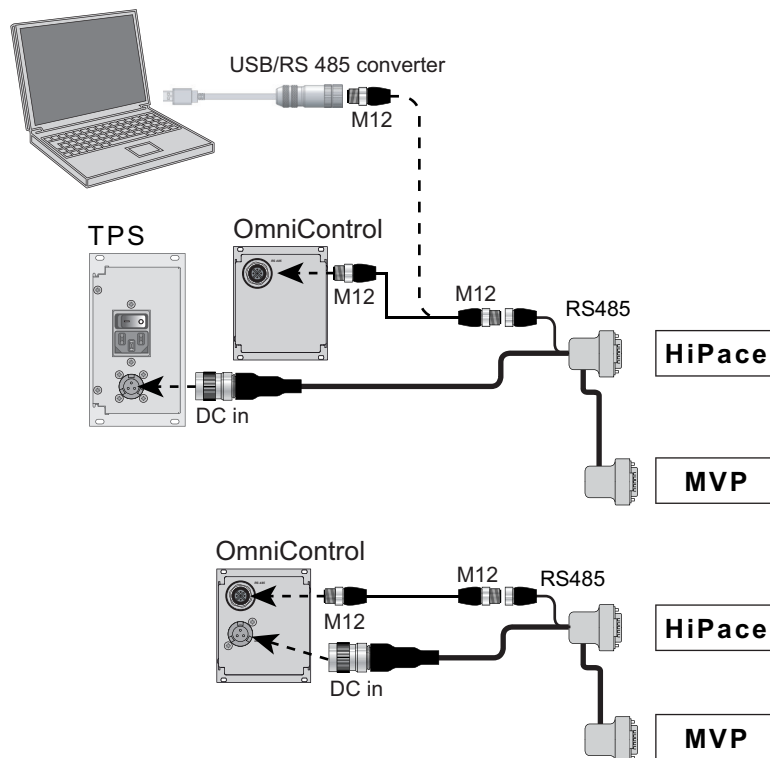


Fig. 8: Connection options via interface RS-485

Connecting control units for Pfeiffer Vacuum or a PC

- ▶ Use the supplied control unit connecting cable or a cable from Pfeiffer Vacuum's accessory range.
- ▶ Connect the control unit to the RS-485 interface.
- ▶ Connect a PC via the USB/RS-485 converter.

6.5 Pfeiffer Vacuum protocol for RS-485 interface

6.5.1 Telegram frame

The telegram frame of the Pfeiffer Vacuum protocol contains only ASCII characters [32; 127], the exception being the end character of the telegram C_R . Basically, a host  (e.g., a PC) sends a telegram, which a unit  (e.g., electronic drive unit or gauge) responds to.

a2	a1	a0	*	0	n2	n1	n0	l1	l0	dn	...	d0	c2	c1	c0	C_R
a2 – a0				Device address 												
				- Individual address of the unit ["001"; "255"] - Group address "9xx" for all identical units (no response) - Global address "000" for all units on the bus (no response)												
*				Action according to telegram description												
n2 – n0				Pfeiffer Vacuum parameter numbers												
l1 – l0				Data length dn to d0												
dn – d0				Data in the respective data type (see chapter "Data types", page 28).												
c2 – c0				Checksum (sum of ASCII values of cells a2 to d0) modulo 256												
C_R				Line break (ASCII 13)												

6.5.2 Telegram description

Data query  -->  ?

a2	a1	a0	0	0	n2	n1	n0	0	2	=	?	c2	c1	c0	C_R
----	----	----	---	---	----	----	----	---	---	---	---	----	----	----	-------

Control command  -->  !

a2	a1	a0	1	0	n2	n1	n0	l1	l0	dn	...	d0	c2	c1	c0	C_R
----	----	----	---	---	----	----	----	----	----	----	-----	----	----	----	----	-------

Data response/control command understood  --> 

a2	a1	a0	1	0	n2	n1	n0	l1	l0	dn	...	d0	c2	c1	c0	C_R
----	----	----	---	---	----	----	----	----	----	----	-----	----	----	----	----	-------

Error message  --> 

a2	a1	a0	1	0	n2	n1	n0	0	6	N	O	_	D	E	F	c2	c1	c0	C_R
										_	Red	A	N	G	E				
										_	L	O	G	I	C				

NO_DEF Parameter number n2 – n0 does not exist
 _RANGE Data dn–d0 outside the permissible range
 _LOGIC Logical access error

6.5.3 Telegram example 1

Data query

Current rotation speed (parameter [P:309], device address: "123")

 -->  ?	1	2	3	0	0	3	0	9	0	2	=	?	1	1	2	C_R
ASCII	49	50	51	48	48	51	48	57	48	50	61	63	49	49	50	13

Data response: 633 Hz

Current rotation speed (parameter [P:309], device address: "123")

 --> 	1	2	3	1	0	3	0	9	0	6	0	0	0	6	3	3	0	3	7	C_R
ASCII	49	50	51	49	48	51	48	57	48	54	48	48	48	54	51	51	48	51	55	13

6.5.4 Telegram example 2

Control command

Switch on the pumping station (parameter **[P:010]**, device address: "042")

 -->  !	0	4	2	1	0	0	1	0	0	6	1	1	1	1	1	1	0	2	0	C _R
ASCII	48	52	50	49	48	48	49	48	48	54	49	49	49	49	49	49	48	50	48	13

Control command understood

Switch on the pumping station (parameter **[P:010]**, device address: "042")

 --> 	0	4	2	1	0	0	1	0	0	6	1	1	1	1	1	1	0	2	0	C _R
ASCII	48	52	50	49	48	48	49	48	48	54	49	49	49	49	49	49	48	50	48	13

6.5.5 Data types

No.	Data type	Description	Length l1 – l0	Example
0	boolean_old	Logical value (false/true)	06	000000 is equivalent to false 111111 is equivalent to true
1	u_integer	Positive whole number	06	000000 to 999999
2	u_real	Fixed point number (unsigned)	06	001571 corresponds with 15.71
4	string	Any character string with 6 characters. ASCII codes between 32 and 127	06	TC_110, TM_700
6	boolean_new	Logical value (false/true)	01	0 is equivalent to false 1 is equivalent to true
7	u_short_int	Positive whole number	03	000 to 999
10	u_expo_new	Positive exponential number. The last of both digits are the exponent with a deduction of 20.	06	100023 is equivalent to $1,0 \cdot 10^3$ 100000 is equivalent to $1,0 \cdot 10^{-20}$
11	string16	Any character string with 16 characters. ASCII codes between 32 and 127	16	BrezelBier&Wurst
12	string8	Any character string with 8 characters. ASCII codes between 32 and 127	08	Example

7 Parameter set

7.1 General

All variables of the vacuum pump relevant for the function are stored as parameters in the electronic drive unit. Each parameter has a three-digit number and a description. Parameters can be used via Pfeiffer Vacuum control units or via RS-485 with the Pfeiffer Vacuum protocol.

Conventions

Parameters are printed in bold as **three-digit numbers** in **square brackets**. The designation can also be specified.

Example: **[P:312]** Software version

Factory settings

The electronic drive unit is pre-programmed in the factory. This makes a more direct and safe operation of the vacuum pump possible without additional configuration.

You can configure the following parameters for the specific process:

- **[P:026]**: rotation speed setting mode on/off.
- **[P:707]**: specification of the rotation speed setting mode with 25 to 134% of the nominal rotation speed.
- **[P:002]**: stand-by mode on/off.

#	Three digit number of the parameter
Indicator	Display of parameter description
Description	Brief description of the parameters
Functions	Function description of the parameters
Data type	Type of formatting of the parameter for the use with the Pfeiffer Vacuum protocol
Access type	R (read): Read access; W (write): Write access
Unit	Physical unit of the described variable
min. / max.	Permissible limit values for the entry of a value
default	Factory default pre-setting (partially pump-specific)
	The parameter can be saved persistently in the electronic drive unit

Tbl. 10: Explanation and meaning of the parameters

7.2 Control commands

#	Indicator	Designations	Functions	Data type	Access type	Unit	min.	max.	de- fault	
002	Stand-by	Stand-by	0 = off 1 = on	0	RW		0	1	0	✓
010	PumpgStatn	Pump	0 = off 1 = on	0	RW		0	1	0	✓
019	Cfg DO2	Output DO2 configuration	1 = No error 2 = Error 5 = Set rotation speed reached 6 = Pump on 9 = "0" 10 = "1" 11 = Remote priority active	7	RW		0	20	6	✓
024	Cfg DO1	Output DO1 configuration	Settings, see [P:019]	7	RW		0	20	1	✓

#	Indicator	Designations	Functions	Data type	Access type	Unit	min.	max.	de-fault	
026	SpdSet-Mode	Rotation speed setting mode	0 = off 1 = on	7	RW		0	1	0	✓
060	CtrlVialnt	Operate via interface	1 = remote 2 = RS-485 4 = PV.can 32 = Keys on the front panel 255 = Interface selection	7	RW		1	255	1	✓
061	IntSelLckd	Interface selection locked	0 = off 1 = on	0	RW		0	1	0	✓

Tbl. 11: Parameter set | Control commands

7.3 Status requests

#	Indicator	Designations	Functions	Data type	Access type	Unit	min.	max.	de-fault	
303	Error code	Error code		4	R					
309	ActualSpd	Actual rotation speed (Hz)		1	R	Hz				
310	DrvCurrent	Drive current		2	R	A	0	9999.99		
311	OpHrsPump	Pump operating hours		1	R	h				✓
312	Fw version	Software version interface circuit board		4	R					
313	DrvVoltage	Supply voltage	Voltage in x.xx V	2	R	V				
315	Nominal Spd	Nominal rotation speed (Hz)		1	R	Hz				
316	DrvPower	Drive power	Output in xW	1	R	W				
330	TempPump	Pump temperature		1	R	°C	0	999999		
349	ElecName	Device name designation		4	R					
352	FWVERSDRV	Firmware of the FC		4	R					
354	HW Version	Hardware version interface circuit board		4	R					
398	ActualSpd	Actual rotation speed (rpm)		1	R	rpm				
399	NominalSpd	Nominal rotation speed (rpm)		1	R	rpm				

Tbl. 12: Parameter set | Status requests

7.4 Reference value inputs

#	Indicator	Designations	Functions	Data type	Access type	Unit	min.	max.	default	
707	SpdSVal	Set value in rotation speed setting mode	Set rotation speed in x.x% of the nominal rotation speed	2	RW	%	25	170	75	
717	StdbySVal	Rotation speed set value in stand-by operation		2	RW	%	30	100	66.7	
797	RS485Adr	RS-485 Interface address		1	RW		1	255	1	

Tbl. 13: Parameter set | Reference value inputs

8 Operation

8.1 Commissioning vacuum pump

WARNING

Danger of poisoning due to toxic process media escaping from the exhaust duct.

During operation with no exhaust duct, the vacuum pump allows exhaust gases and vapors to escape freely into the air. There is a risk of injury and fatalities due to poisoning in processes with toxic process media.

- ▶ Observe the pertinent regulations for handling toxic process media.
- ▶ Safely purge toxic process media via an exhaust duct.
- ▶ Use appropriate filter equipment to separate toxic process media.

NOTICE

Vacuum pump damage caused by overpressure

Mixing up the connections causes overpressure overload. The vacuum pump can be started only against max. 1 bar differential pressure between inlet and outlet; otherwise the motor jams and sustains damage.

- ▶ Make sure that the distributor is installed correctly on the diaphragm heads, before connecting the vacuum pump to the vacuum equipment.
- ▶ Before commissioning, make completely sure that the pressure side pressure is below the permissible limit.

Parameter	Name	Designation	Adjustment, setting
[P:707]	SpdSVal	Speed-control operation specification	75 %

Tbl. 14: Factory setting of the electronic drive unit for diaphragm pumps when delivered

Procedure

1. Compare the frequency specifications on the rating plate with the available supply voltage.
2. Check the exhaust connection for free passage (max. permissible pressure: 1100 hPa absolute).
3. Actuate the shut-off units so that the shut-off units open before or at the same time as the vacuum pump starts up.

8.2 Switching on the vacuum pump

When pumping down dry gases, no special precautions are required.

CAUTION

Danger of burns on hot surfaces

Depending on the operating and ambient conditions, the surface temperature of the vacuum pump can increase to above 70 °C.

- ▶ Provide suitable touch protection.



Warm-up time of the vacuum pump

The warm-up time depends on the ambient temperature and takes at least 30 minutes.

Prerequisite

- required cable connections established

Switching on vacuum pump

1. If required, switch the vacuum pump on in each pressure range in one of the 3 possible ways:
 - using a mains switch on the power supply pack with respective bridges on the 15-pin D-Sub socket
 - via a control unit, e.g. OmniControl or PC
 - using the PLC control unit on the 15-pin D-Sub socket
2. Allow the vacuum pump to warm up prior to process start, with the vacuum connection closed.

The vacuum pump achieves the specified throughput and final pressure values once the operating temperature has been reached.

8.3 Configuring connections with Pfeiffer Vacuum parameter set

The electronic drive unit is pre-configured with the factory default basic functions and is ready for operation. For individual requirements, most connections for the electronic drive unit can be configured with the parameter set.

8.3.1 Configuring the digital outputs

Option	Description
1 = No error	active, with trouble-free operation
2 = Error	active, if the error message is active
5 = Set rotation speed reached	active, once the set rotation speed is reached
6 = Pump on	active, if pump on, motor on and no error
9 = always 0	GND for the control of an external device
10 = always 1	V+ for the control of an external device
11 = Remote	active, if the remote priority is active

Tbl. 15: Configure parameters [P:019] and [P:024]

Procedure

- Perform the configuration according to the table.

8.3.2 Selecting the interfaces

Option [P:060]	Description
1 = remote	Operation via connection "remote"
2 = RS-485	Operation via connection "RS-485"
4 = PV.can	For service purposes only
32 = Keys on the front panel	Manual operation
Option [P:061]	
0 = off	Interface selection can be set via [P:060].
1 = on	Interface selection locked

Tbl. 16: Parameters [P:060] and [P:061]

Procedure

- Select the interface according to the table.

8.4 Operating modes

- Operating without control unit
- Operation via an external control unit
- Operation via RS-485 and Pfeiffer Vacuum control unit or PC

The connection of a Pfeiffer Vacuum control unit permits the controlling of the vacuum pump via the parameters fixed in the electronic drive unit.

**Automatic start**

After bypassing the contacts at pins 2, 7, 5 on the connecting plug or when using a connection cable with respective bridges and applying the supply voltage, the vacuum pump starts up immediately. Pfeiffer Vacuum recommends switching on the voltage supply immediately before operation.

Operating vacuum pump without controller

1. Connect the respective connecting cable with bridges to the "remote" connection of the electronic drive unit.
2. Provide the current supply for the power supply pack.
3. Switch on the current supply on the power supply pack.

After applying the operating voltage, the vacuum pump carries out a self-test to check the supply voltage. After a successfully completed self-test, the vacuum pump starts.

Operating vacuum pump with external controller

1. Connect a remote control via the "remote" 15-pin D-Sub socket.
2. Control the vacuum pump by means of the "PLC level".

Operating vacuum pump with control unit

1. Observe the operating instructions of the control unit.
2. Connect the respective control unit at the "remote" connection.
3. Switch on the current supply on the power supply pack or on the control unit.
4. Make the desired settings via the RS-485 using the control unit.

8.4.1 Rotation speed setting operation

The set rotation speed is selected via the parameter **[P:707]** in the range **30 to 134%** of the set rotation speed. If the speed is higher than the set rotation speed (> 100%), an increased pump wear is possible.

The rotation speed setting mode has priority over the stand-by mode.

**Permissible rotation speed range of the vacuum pump**

Parameterization in rotation speed setting mode is subject to the permissible rotation speed range of the respective vacuum pump. The electronic drive unit regulates automatically to the next valid value.

Set the rotation speed setting mode

1. Set the parameter **[P:707]** to the required value in %.
2. Set the parameter **[P:026]** to "1".
3. Check the set rotation speed (parameter **[P:308]** or **[P:397]**).

8.4.2 Standby mode

Pfeiffer Vacuum recommends standby mode for during process and production stops. With active stand-by mode, the electronic drive unit reduces the speed of the vacuum pump in the range of **30 to 100%** of the nominal speed. The factory setting is 66.7% of the nominal speed.

Setting the related parameters

1. Set the parameter **[P:717]** to the required value in %.
2. Set the parameter **[P:026]** to "0".
3. Set the parameter **[P:002]** to "1".
4. Check the set rotation speed (parameter **[P:308]** or **[P:397]**).

8.4.3 Normal operation

The vacuum pump starts in boost mode for max. 5 minutes with the maximum speed. The current is monitored at the same time. If the current is below a threshold value, the speed is controlled down to nominal rotation speed. In case the power input increases again, automatic increase of the speed does not take place. An increase in speed is reached again after vacuum pump stop / start.

Setting the related parameters

1. Set the parameter **[P:002]** to "0".
2. Set the parameter **[P:026]** to "0".
3. Check the set rotation speed (parameter **[P:308]**).

8.5 Pumping condensable vapors

Vapors or moisture from pumped media impair the throughput after condensation in the vacuum pump. Using the optional flushing gas nozzle improves the discharge of condensate, and the pump achieves the specified final vacuum more quickly.

⚠ WARNING

Risk of explosion from reactive, explosive or other hazardous gas / air mixtures.

Uncontrolled inlet of gas at the flushing gas nozzle leads to sparks generated mechanically following diaphragm rupture, hot surfaces or to explosive gas / air mixtures in the vacuum system caused by static electricity.

- If necessary, use inert gas for supplying the flushing gas in order to avoid a potential ignition.

⚠ CAUTION

Risk of poisoning due to toxic process gases escaping

For intake pressures > 500 hPa, process gas may escape from the flushing gas nozzle.

- If necessary, use an extraction line.



Maximum inlet pressure

The maximum permissible inlet pressure at the flushing gas nozzle is 1,200 hPa absolute.

Accessories required

- Flushing gas nozzle with filter (order number PK 050 304 -U)

Required tools

- Allen key, **WAF 5**
- Open-end wrench, **WAF 14**

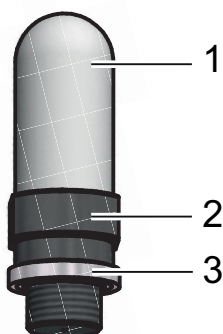


Fig. 9: Flushing gas nozzle with filter

- | | |
|--|-------------------|
| 1 Filter (1/8" thread) | 3 VS sealing ring |
| 2 Nozzle adapter for inert gas connection (1/8" thread, nozzle Ø 0.2 mm) | |

Mounting the flushing gas nozzle

1. Unscrew and remove the sealing screw from the vacuum pump.
2. Screw the flushing gas nozzle into the vacuum pump.
3. Manually unscrew and remove the filter from the nozzle adapter to connect an inert gas line.

8.6 Operation monitoring

8.6.1 Operating mode display via LED

The LED on the electronic drive unit indicates the basic operating conditions. A differentiated error and warning display is only possible for operation with a Pfeiffer Vacuum control unit.

Indicator	Activity	Meaning
	None	No adequate power supply
	10% active, 1 Hz	- No error - Vacuum pump "OFF" - Vacuum pump stationary
	50% active, 1 Hz	- No error - Vacuum pump "OFF" - Vacuum pump still rotating
	90% active, 1 Hz	- No error - Vacuum pump "ON" - Set rotation speed not reached
	100% active	- No error - Vacuum pump "ON" - Set rotation speed reached
	50% active, 10, Hz	Error

Tbl. 17: Behavior and meaning of the LED of the electronic drive unit

8.6.2 Temperature monitoring

If the pump temperature becomes too high ($> 75\text{ °C}$), the electronic drive unit reduces the motor rotation speed to the nominal rotation speed to prevent the vacuum pump from overheating.

After cooling down ($< 72\text{ °C}$), the vacuum pump starts to run at set rotation speed.

8.7 Switching off the vacuum pump

Procedure

1. Allow the vacuum pump to run on for 5 to 10 minutes with the vacuum connection open to allow any condensate that may be present to be removed from the vacuum pump.
2. If required, switch the vacuum pump off in each pressure range.
3. Disconnect the voltage supply on the power supply pack.

9 Maintenance

9.1 Maintenance instructions

WARNING

Health hazard through poisoning from toxic contaminated components or devices

Toxic process media result in contamination of devices or parts of them. During maintenance work, there is a risk to health from contact with these poisonous substances. Illegal disposal of toxic substances causes environmental damage.

- ▶ Take suitable safety precautions and prevent health hazards or environmental pollution by toxic process media.
- ▶ Decontaminate affected parts before carrying out maintenance work.
- ▶ Wear protective equipment.

CAUTION

Danger of injury from moving parts

After a power failure or a standstill as a result of overheating, the motor restarts automatically. There is a risk of injury to fingers and hands if they enter the operating range of rotating parts.

- ▶ Safely disconnect motor from the mains.
- ▶ Secure the motor against reactivation.
- ▶ Dismantle the vacuum pump for inspection, away from the system if necessary.

NOTICE

Damage caused by unsuitable cleaning agents

Unsuitable cleaning agents damage pump parts.

- ▶ Use only approved cleaning agents to clean pump parts.
- ▶ Use only a clean dry cloth to clean the diaphragms and valves.
- ▶ Do not use any alcohol or other cleaning agents to clean the diaphragms and valves.

NOTICE

Danger of property damage from improper maintenance

Unprofessional work on the vacuum pump will lead to damage for which Pfeiffer Vacuum accepts no liability.

- ▶ We recommend taking advantage of our service training program.
- ▶ When ordering spare parts, specify the information on the nameplate.

The following section describes the tasks for cleaning and maintaining the vacuum pump.

Valves and diaphragms are wear parts.

Component	Operating hours
Diaphragms	15 000
Valves	15 000
Motor bearing	40 000

Tbl. 18: Typical service life with normal use

Prerequisites

- Vacuum pump switched off
- Vacuum pump vented to atmospheric pressure
- Vacuum pump cooled

Required materials

- Cloth (clean, lint-free)
- Water or mild soap solution

Servicing the vacuum pump

1. Disconnect the vacuum pump from the voltage supply.
2. Secure the motor against reactivation.
3. Remove any external contamination on the vacuum pump with cloth slightly moistened with water or a mild soap solution.
4. Allow all cleaned parts to dry well.
5. For maintenance work, only dismantle the vacuum pump to the extend needed.
6. Clean the suction chamber, diaphragms, and valves with a dry cloth.
7. Examine the diaphragms and valves for cracks at the latest when the pressure values achieved decrease.

9.2 Checklist for inspection and maintenance**Maintenance frequency and service lives**

Maintenance frequency and service lives are process-dependent. Chemical and thermal loads, as well as contamination, shorten the recommended reference values.

- Determine the specific service lives during the first operating interval.
- Please contact Pfeiffer Vacuum Customer Service if you wish to shorten the maintenance intervals.

**Maintenance by Pfeiffer Vacuum Service**

We recommend that Pfeiffer Vacuum Service carry out maintenance work. If the specified intervals are exceeded, or if maintenance work is carried out improperly, no warranty or liability claims are accepted on the part of Pfeiffer Vacuum. This also applies wherever parts other than original spare parts are used.

Action	Inspection	Maintenance	Required material
Interval	as required; at least once every six months	as required; at least every 2 years	
Check silencer for contamination ²⁾	■		
Test vacuum pump optically and acoustically	■		
Read out and analyze pump data ³⁾		■	
Clean the vacuum pump		■	
Change the diaphragms and valves		■	Overhaul kit
Change silencer ⁴⁾		■	Silencer
Performing a function test		■	
Performing incoming inspection		■	
Clean vacuum pump completely		■	
Change gas ballast filter ⁵⁾		■	Filter
Replace wear parts		■	Diaphragms, valves, sealing rings, silencers

Tbl. 19: Maintenance intervals

9.3 Dismantling / installing the handle**Required tools**

- Phillips head (crosshead) screwdriver, **size 2**

2) if present

3) DC pumps only

4) if present

5) if present

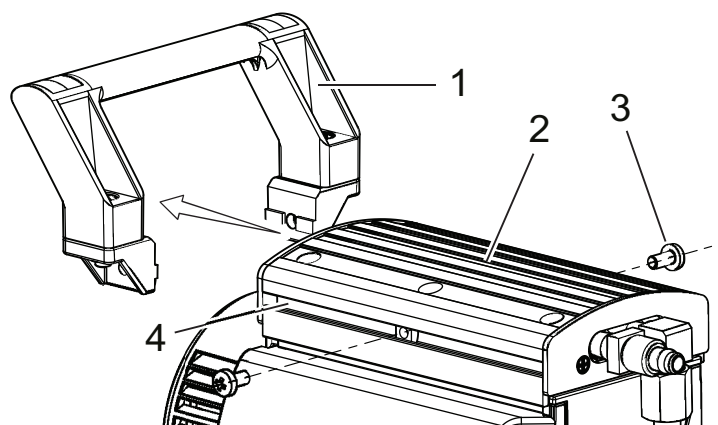


Fig. 10: Handle

- | | |
|-----------------|--------------|
| 1 Handle | 3 Screw (2×) |
| 2 Housing cover | 4 Groove |

Dismantling the handle

1. Loosen the screws of the handle.
2. Push the handle from the groove in the housing cover.

Installing the handle

1. Push the handle into the groove in the housing cover.
2. Pay attention to the correct position of the handle above the center of gravity of the vacuum pump.
3. Fasten the screws of the handle to the nuts in the housing cover.

9.4 Change the diaphragms and valves

NOTICE

Property damage from incorrect installation

Change in dead volume due to incorrect installation of the original spacer disks impairs the final vacuum or leads to bearing damage.

- During disassembly, keep the existing spacer disks separate per installation location.
- Reinstall the same number of original spacer disks per diaphragm head.

9.4.1 Dismantle the pump head and valves



Avoid interchanging components

Always dismantle and mount only the pump head on one pump side in order to avoid interchanging the components.

The procedures are identical for both pump heads.

Prerequisites

- Handle dismantled
- Diaphragm pump supported

Required tools

- Open-end wrench, **WAF 15**
- Open-end wrench, **WAF 16**
- Allen key, **WAF 4**

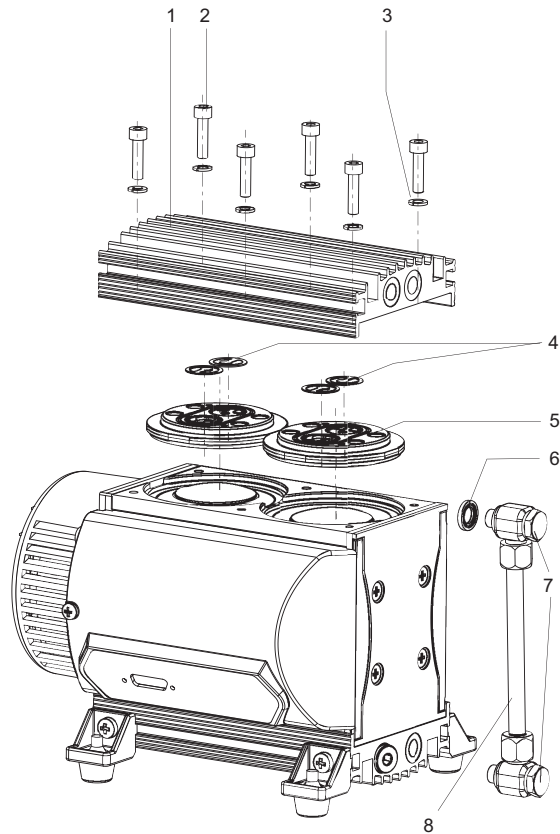


Fig. 11: Pump head and valves

- | | |
|-----------------------|---------------------|
| 1 Housing cover 1 | 5 Head cover (2×) |
| 2 Cylinder screw (6×) | 6 Sealing ring (2×) |
| 3 Lock washer (6×) | 7 Banjo bolt (2×) |
| 4 Valve (4×) | 8 Hose connection |

Procedure

1. Loosen the screw fitting.
2. Turn the elbow fitting max. 1/4 revolution anti-clockwise, until the hose connection can be loosened.
3. Loosen the screw fitting.
4. Dismantle the hose connection.
5. Unscrew the cylinder screws on the housing cover.
6. Remove the housing cover.
7. Remove the head cover each with both valves.
 - If the valves stick to the housing cover, carefully release the valves; otherwise, remove the valves from the head covers.

9.4.2 Dismantling the diaphragms

Prerequisite

- Pump head and valves dismantled

Required tool

- Diaphragm key **WAF 46**

Spare parts required

- Inspection set, including diaphragm key

Required consumable materials

- Clean, dry cloth
- Alcohol or cleaning solvent, if necessary

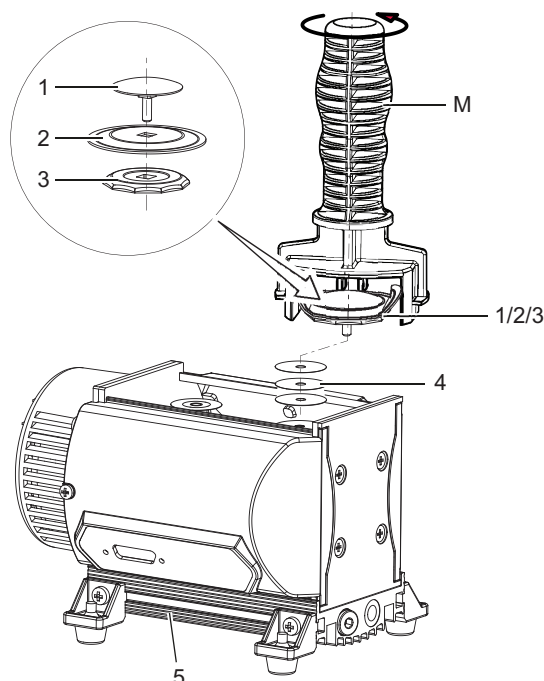


Fig. 12: Replacing the diaphragms

- | | |
|---------------------|-----------------|
| 1 Clamping washer | 4 Spacer disks |
| 2 Diaphragm | 5 Pump head 2 |
| 3 Supporting washer | M Diaphragm key |

Dismantling the diaphragms

- Carefully bend the diaphragm up on the side and while it is bent, attach the diaphragm key to the supporting washer.
 - Do not use a sharp-edged tool.
- Unscrew the supporting washer, diaphragm, and clamping washer.
- Pay attention to the spacer disks under the diaphragm.
 - Reuse the original spacer disks for each installation location when remounting.
- Remove the supporting washer and the diaphragm from the square of the clamping washer connecting screw.
 - If the diaphragm is stuck to the supporting washer, use alcohol or cleaning solvent to release it.

9.4.3 Installing the diaphragm

Required tools

- Diaphragm key **WAF 46**

Spare parts required

- Overhaul kit incl. diaphragm key

Required consumables

- Clean, dry cloth
- Alcohol or cleaning solvent, if necessary

Installing the diaphragm

- Clean all parts with a dry, clean cloth.
- Replace the wearing parts (valves and diaphragms) as per the spare part package.
- Place the diaphragm (blue side up) and the supporting washer on the square of the clamping washer connecting screw.
 - Pay attention to the seating of the diaphragm in the square of the connecting screw.
- Place the original spacer disks on the respective connection rod.
- Tighten the supporting washer with the diaphragm key on the connection rod.

9.4.4 Installing the pump head and valves



Avoid interchanging components

Always dismantle and mount only the pump head on one pump side in order to avoid interchanging the components.

The procedures are identical for both pump heads.

Prerequisite

- Diaphragms installed

Required tools

- Open-end wrench, **WAF 15**
- Open-end wrench, **WAF 16**
- Allen key, **WAF 4**
- Calibrated torque wrench (tightening factor ≤ 1.6)

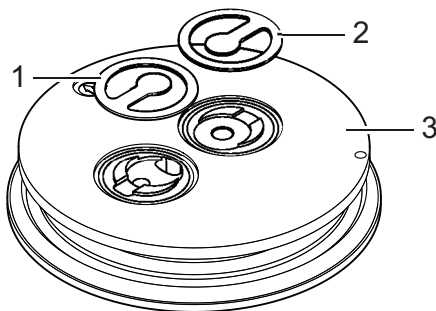


Fig. 13: Correct mounting orientation of the valves in the head cover

- | | |
|------------------|--------------|
| 1 Valve (inlet) | 3 Head cover |
| 2 Valve (outlet) | |

Procedure

1. Insert the valves into the head cover.
 - Make sure that the valves have the correct mounting orientation.
2. Mount the housing cover.
3. First, screw on the cylinder screws of the housing cover loosely by hand, alternating screws in a diagonal pattern.
4. Then tighten the cylinder screws of the housing cover, alternating screws in a diagonal pattern.
 - Tightening torque: **max. 6.0 ± 1 Nm**
5. Mount the hose connection and the banjo bolts.

9.5 Checking final pressure

The leakage rate of the vacuum pump ensures that the specified final pressure is reached and prevents the formation of potentially explosive mixtures in the interior of the vacuum pump

After the diaphragm or valve is replaced, the vacuum pump reaches the specified final pressure value only after a run-in period of several hours.



Test values

Observe the separate test specification.

- with gas ballast: < 3 hPa
- without gas ballast: < 2.0 hPa

Procedure

1. Measure the final pressure of the vacuum pump.
 - after work on the system, e.g., maintenance
 - before the vacuum pump is reinserted
2. Compare the measured final pressure with the test values and the specifications in the technical data.

Prerequisites

- Vacuum pump mounted
- Vacuum pump correctly installed electrically

Required aids

- Vacuum chamber (3 liter)
- Pressure gauge
- Flushing gas nozzle (optional)

Measuring final pressure

1. Connect a vacuum chamber (volume approx. 3 l) to a vacuum connection.
2. Also connect a pressure gauge (e.g. Pirani gauge) to the vacuum chamber on the vacuum side.
3. Operate the vacuum pump with open flushing gas nozzle (if available) for at least 2 hours.
4. If unusual noises develop, immediately switch off the vacuum pump and check the position of the clamping washers.
5. Measure the final pressure.

Comparing the ultimate pressure with the test value

1. Compare the measured ultimate pressure with the test value.
2. If the values differ significantly from the specified values after replacing the diaphragm and valve, and if there are no changes after a run-in period, perform the following steps:
 - Check the compression couplings of the connection hoses to the pump heads.
 - Check the valve seats and the suction chambers.
 - Repeat the ultimate pressure measurement.

10 Decommissioning

Before shutting down the vacuum pump, observe the following instructions to adequately protect the interior of the vacuum pump (suction chamber) from corrosion:

Procedure for temporary vacuum pump shutdowns

1. Allow the vacuum pump to run on for 5 to 10 minutes with the vacuum connection open to allow any condensate that may be present to be removed from the vacuum pump.
2. If media has entered the vacuum pump that may attack the pump materials or lead to deposits, clean the inside of the pump heads.

Procedure for longer vacuum pump shutdowns

1. Disconnect the vacuum pump from the vacuum system.
2. Remove the vacuum pump from the vacuum system if necessary.
3. Seal the vacuum connection with the original sealing plugs.
4. Store the vacuum pump in dry, dust-free rooms, within the specified ambient conditions.
5. Pack the vacuum pump together with a drying agent in a plastic bag, and seal the vacuum pump airtight if it is to be stored in rooms with damp or aggressive atmospheres.

11 Recycling and disposal

⚠ WARNING

Health hazard through poisoning from toxic contaminated components or devices

Toxic process media result in contamination of devices or parts of them. During maintenance work, there is a risk to health from contact with these poisonous substances. Illegal disposal of toxic substances causes environmental damage.

- ▶ Take suitable safety precautions and prevent health hazards or environmental pollution by toxic process media.
- ▶ Decontaminate affected parts before carrying out maintenance work.
- ▶ Wear protective equipment.



Environmental protection

You **must** dispose of the product and its components in accordance with all applicable regulations for protecting people, the environment and nature.

- Help to reduce the wastage of natural resources.
- Prevent contamination.

11.1 General disposal information

Pfeiffer Vacuum products contain materials that you must recycle.

- ▶ Dispose of our products according to the following:
 - Iron
 - Aluminium
 - Copper
 - Synthetic
 - Electronic components
 - Oil and fat, solvent-free
- ▶ Observe the special precautionary measures when disposing of:
 - Fluoroelastomers (FKM)
 - Potentially contaminated components that come into contact with media

11.2 Dispose of diaphragm pumps

Pfeiffer Vacuum diaphragm pumps contain materials that you must recycle.

1. Disconnect the electronic drive unit.
2. Dismantle the motor.
3. Decontaminate the components that come into contact with process gases.
4. Separate the components into recyclable materials.
5. Recycle the non-contaminated components.
6. Dispose of the product or components in a safe manner according to locally applicable regulations.

12 Malfunctions

12.1 General

CAUTION

Danger of injury from moving parts

After a power failure or a standstill as a result of overheating, the motor restarts automatically. There is a risk of injury to fingers and hands if they enter the operating range of rotating parts.

- ▶ Safely disconnect motor from the mains.
- ▶ Secure the motor against reactivation.
- ▶ Dismantle the vacuum pump for inspection, away from the system if necessary.

CAUTION

Danger of burns on hot surfaces

In the event of a fault, the surface temperature of the vacuum pump can increase to above 105°C.

- ▶ Allow the vacuum pump to cool down before carrying out any work.
- ▶ Wear personal protective equipment if necessary.

NOTICE

Danger of property damage from improper maintenance

Unprofessional work on the vacuum pump will lead to damage for which Pfeiffer Vacuum accepts no liability.

- ▶ We recommend taking advantage of our service training program.
- ▶ When ordering spare parts, specify the information on the nameplate.

Vacuum pump and electronic drive unit malfunctions always result in a warning or error message. In both cases, you will receive an error code that you can read out via the interface of the electronic drive unit. If an error occurs, the vacuum pump switches off.

Should malfunctions occur, you can find information about potential causes and how to fix them here:

Problem	Possible causes	Remedy
Vacuum pump will not start up	No supply voltage or voltage does not correspond to the motor data	Check the supply voltage.
	Pump temperature too low	Heat the vacuum pump to a temperature of > 5°C.
	Thermal protection of the motor has triggered	Identify and eliminate the cause and allow vacuum pump to cool if necessary.
	Diaphragms or valves dirty	Clean the vacuum pump.
	Overpressure in the exhaust lead	If required, check and clean the exhaust line.
Vacuum pump switches off after a while after being started	Thermal protection of the motor has triggered	Identify and eliminate the cause and allow vacuum pump to cool if necessary.
	Exhaust pressure too high	Check exhaust line outlet opening and exhaust side accessories.

Problem	Possible causes	Remedy
Vacuum pump does not reach the specified ultimate pressure	Condensate in the vacuum pump	Operate the vacuum pump at atmospheric pressure for a longer period.
	Flushing gas nozzle/gas ballast valve open	Close the flushing gas nozzle/the gas ballast valve.
	Soiled valves or diaphragms	If necessary, clean or replace the valves and diaphragms.
	Leak in system	Locate and eliminate the leak.
Pumping speed of vacuum pump too low	The intake line is not suitably dimensioned	Make sure that connections are short and cross sections adequately dimensioned.
	Exhaust pressure too high	Check exhaust line outlet opening and exhaust side accessories.
Unusual noises during operation	Defective valves or diaphragms	If necessary, clean or replace the valves and diaphragms.
	Soiled suction chamber	Clean suction chamber.
	The silencer is loose or missing	If necessary, check or replace the silencer.
	Valves dirty or defective	If necessary, clean or replace the valves and diaphragms.
	Connection rod or motor bearing defective	Contact Pfeiffer Vacuum Service.

Tbl. 20: Troubleshooting on diaphragm pumps

12.2 Error codes

Errors (** Error E—— **) always cause the vacuum pump to be switched off.

Warnings (* Warning F —— *) do not cause the vacuum pump to be switched off.

Handling malfunction messages

1. Read out error codes via Pfeiffer Vacuum control units or a PC.
2. Remove the cause of the malfunction.
3. Reset the malfunction message with parameter **[P:009]**.
 - Use preconfigured interfaces or screen tiles on Pfeiffer Vacuum control units.

Error code	Problem	Possible causes	Remedy
Err098	Internal communication error between interface board and drive	–	Contact Pfeiffer Vacuum Service.
Err117	Vacuum pump excess temperature (> 90 °C)	Insufficient cooling	- Improve the cooling - Check the operating conditions
Err174	Vacuum pump blocked	–	Contact Pfeiffer Vacuum Service.

Tbl. 21: Error messages for vacuum pump

Error code	Problem	Possible causes	Remedy
Wrn117	High temperature vacuum pump (> 75 °C)	Insufficient cooling	- Improve the cooling - Check the operating conditions

Tbl. 22: Warning messages for vacuum pump

13 Service solutions by Pfeiffer Vacuum

We offer first-class service

High vacuum component service life, in combination with low downtime, are clear expectations that you place on us. We meet your needs with efficient products and outstanding service.

We are always focused on perfecting our core competence – servicing of vacuum components. Once you have purchased a product from Pfeiffer Vacuum, our service is far from over. This is often exactly where service begins. Obviously, in proven Pfeiffer Vacuum quality.

Our professional sales and service employees are available to provide you with reliable assistance, worldwide. Pfeiffer Vacuum offers an entire range of services, from [original replacement parts](#) to [service contracts](#).

Make use of Pfeiffer Vacuum service

Whether preventive, on-site service carried out by our field service, fast replacement with mint condition replacement products, or repair carried out in a [Service Center](#) near you – you have various options for maintaining your equipment availability. You can find more detailed information and addresses on our homepage, in the section.

You can obtain advice on the optimal solution for you, from your [Pfeiffer Vacuum representative](#).

For fast and smooth service process handling, we recommend the following:



1. Download the up-to-date form templates.
 - [Explanations of service requests](#)
 - [Service requests](#)
 - [Contamination declaration](#)



- a) Remove and store all accessories (all external parts, such as valves, protective screens, etc.).
- b) If necessary, drain operating fluid/lubricant.
- c) If necessary, drain coolant.
2. Complete the service request and contamination declaration.



3. Send the forms by email, fax, or post to your local [Service Center](#).




4. You will receive an acknowledgment from Pfeiffer Vacuum.


PFEIFFER VACUUM

Submission of contaminated products


No microbiological, explosive, or radiologically contaminated products will be accepted. Where products are contaminated, or the contamination declaration is missing, Pfeiffer Vacuum will contact you before starting service work. Depending on the product and degree of pollution, **additional decontamination costs** may be incurred.

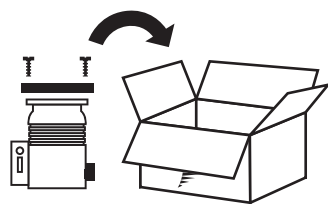

5. Prepare the product for transport in accordance with the provisions in the contamination declaration.
 - a) Neutralize the product with nitrogen or dry air.
 - b) Seal all openings with blind flanges, so that they are airtight.
 - c) Shrink-wrap the product in suitable protective foil.
 - d) Package the product in suitable, stable transport containers only.
 - e) Maintain applicable transport conditions.

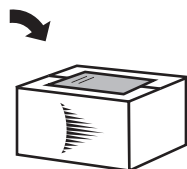

6. Attach the contamination declaration to the **outside** of the packaging.

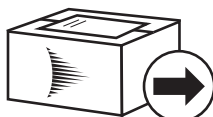

7. Now send your product to your local Service Center.



8. You will receive an acknowledgment/quotation, from Pfeiffer Vacuum.

PFEIFFER VACUUM

Our sales and delivery conditions and repair and maintenance conditions for vacuum devices and components apply to all service orders.

14 Spare part packages

Ordering spare part packages

- ▶ Have the vacuum pump part number to hand, along with other details from the rating plate if necessary.
- ▶ Install original spare parts only.
- ▶ When ordering the inspection set, observe the respective part number of the diaphragm pump.

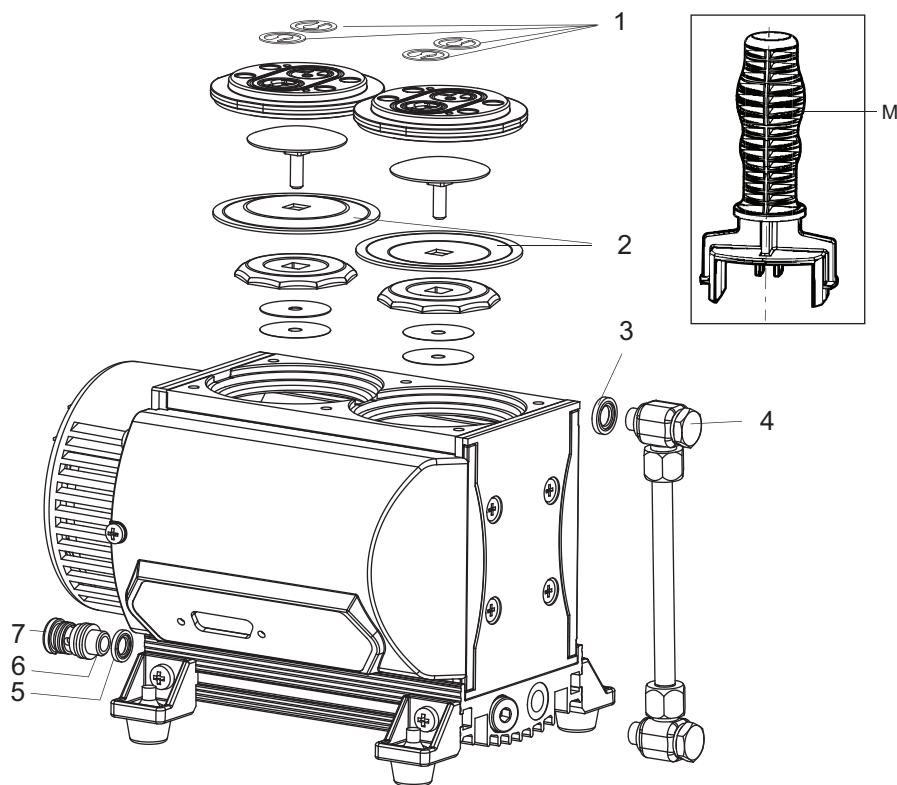


Fig. 14: Spare parts

- | | |
|----------------|-----------------|
| 1 Valves | 5 Seal |
| 2 Diaphragm | 6 Silencer |
| 3 Sealing ring | 7 Hose piece |
| 4 Banjo bolt | M Diaphragm key |

Spare part package	Order number	comprising the following parts
Inspection set, including diaphragm key	PU E22 003 -T for part numbers PK T01 500 and PK T01 501	Pos. 1 (8×) Pos. 2 (4×) Pos. 7 (1×) Pos. M
	PU E22 034 -T for part number PK T01 502	Pos. 1 (8×) Pos. 2 (4×) Pos. M
Sealing ring	P 3529 133 -P	Pos. 3
Banjo bolt	P 0995 943	Pos. 4
Silencer incl. hose piece	P 0995 942	Pos. 6, 7
Diaphragm key	P 0995 941	Pos. M

Tbl. 23: Spare part packages

Spare part package	Order number	comprising the following parts
Filter	P 0105 405	Flushing gas nozzle filter

Tbl. 24: Spart part package flushing gas nozzle

15 Accessories

15.1 Accessory information

Control units

Control units are used to check and adjust operating parameters.

Power supply packs

Power supply packs for fastening to the wall and standard rails or for rack assembly serve the voltage supply.

Cable and adapter

Mains, interface, connection, and extension cables provide a secure and suitable connection. Different lengths on request

Hose connections

Hose connections with adapters for connecting to a turbopump

Screw-in flange

Screw-in flange, including seal for the intake side of the vacuum pump

Flushing gas connection

Avoids moisture accumulating in the vacuum pump

15.2 Ordering accessories



OmniControl variants

You can find additional OmniControl variants on the [Pfeiffer Vacuum website](#).

Selection field	Part number
OmniControl 200, rack unit with integrated power supply pack , No Data, No Gauge/IO Option	PE D50 000 0
OmniControl 200, table unit with integrated power supply pack , No Data, No Gauge/IO Option	PE E50 000 0
TPS 180, power supply pack for wall/standard rail installation	PM 061 341 -T
Interface cable, M12 m straight / M12 m straight, 3 m	PM 061 283 -T
Connecting cable from 24V/48V power supply pack to electronic drive unit. With RS-485 interface	PM 061 350 -T
Connection cable MVP-TC-TPS, 3 m	PE 100 013 -T
USB RS-485 converter	PM 061 207 -T
Y-connector M12 for RS-485	P 4723 010
Hose connection DN 6 x 400 mm with straight connector 1/8" thread and 1/4" thread including seal	P 0920 739
Hose connection DN 6 x 1000 mm with straight connector 1/8" thread and 1/4" thread including seal	P 0920 817
Hose DN 6; (ø 8/6 mm) polyethylene	P 0991 939
Push-in fitting G 1/8" including seal for hose connection (8/6 mm)	P 4131 029
Screwing flange DN 16 ISO-KF / 1/8" thread incl. seal	PK 050 108 -T
Flushing gas nozzle with filter for MVP 015-4 / MVP 020-3 / MVP 030-3	PK 050 304 -U

Tbl. 25: Accessories

16 Technical data and dimensions

16.1 General

Basis for the technical data of Pfeiffer Vacuum diaphragm pumps:

- Specifications according to PNEUROP committee PN5
- ISO 21360:2012: "Vacuum technology - Standard methods for measuring vacuum-pump performance - General description"

The following harmonized standards are fulfilled:

- IEC 61010-1
- UL 61010-1
- CSA 61010-1

	mbar	bar	Pa	hPa	kPa	Torr mm Hg
mbar	1	$1 \cdot 10^{-3}$	100	1	0.1	0.75
bar	1000	1	$1 \cdot 10^5$	1000	100	750
Pa	0.01	$1 \cdot 10^{-5}$	1	0.01	$1 \cdot 10^{-3}$	$7.5 \cdot 10^{-3}$
hPa	1	$1 \cdot 10^{-3}$	100	1	0.1	0.75
kPa	10	0.01	1000	10	1	7.5
Torr mm Hg	1.33	$1.33 \cdot 10^{-3}$	133.32	1.33	0.133	1

1 Pa = 1 N/m²

Tbl. 26: Conversion table: Pressure units

	mbar l/s	Pa m ³ /s	sccm	Torr l/s	atm cm ³ /s
mbar l/s	1	0.1	59.2	0.75	0.987
Pa m ³ /s	10	1	592	7.5	9.87
sccm	$1.69 \cdot 10^{-2}$	$1.69 \cdot 10^{-3}$	1	$1.27 \cdot 10^{-2}$	$1.67 \cdot 10^{-2}$
Torr l/s	1.33	0.133	78.9	1	1.32
atm cm ³ /s	1.01	0.101	59.8	0.76	1

Tbl. 27: Conversion table: Units for gas throughput



Performance data

- The vacuum pump delivers its highest pumping speed at maximum rotation speed.
- The vacuum pump reaches its lowest final pressure in a process-specific range below the nominal rotation speed.

16.2 Technical data

Type designation	MVP 030-3 DC	MVP 030-3 DC	MVP 030-3 DC	MVP 030-3 DC
Part number	PK T01 500	PK T01 501	PK T01 502	PK T01 503
Connection flange (in)	G 1/8"	G 1/8" thread FESTO push-in fitting NPQH-L- G18-Q8-P10	G 1/8"	G 1/8"
Connection flange (out)	G 1/8" with silencer	G 1/8" with silencer	G 1/8"	G 1/8" with silencer
Pumping speed	1.8 m ³ /h	1.8 m ³ /h	1.8 m ³ /h	1.8 m ³ /h
Ultimate pressure	2 hPa	2 hPa	2 hPa	2 hPa
Ultimate pressure with gas ballast	3 hPa	3 hPa	3 hPa	3 hPa
Intake pressure max.	1100 hPa	1100 hPa	1100 hPa	1100 hPa

Type designation	MVP 030-3 DC	MVP 030-3 DC	MVP 030-3 DC	MVP 030-3 DC
Exhaust pressure, max.	1100 hPa	1100 hPa	1100 hPa	1100 hPa
Integral leak rate	$1 \cdot 10^{-2}$ Pa m ³ /s	$1 \cdot 10^{-2}$ Pa m ³ /s	$1 \cdot 10^{-2}$ Pa m ³ /s	$1 \cdot 10^{-2}$ Pa m ³ /s
Operating voltage: DC	24 V	24 V	24 V	24 V
Input voltage: tolerance	±10 %	±10 %	±10 %	±10 %
Rated current consumption	1.5 A	1.5 A	1.5 A	1.5 A
Current, max.	7 A	7 A	7 A	6 A
Rotation speed	600 – 2400 rpm	600 – 2400 rpm	600 – 2400 rpm	600 – 2400 rpm
I/O interfaces	RS-485	RS-485	RS-485	RS-485
Cooling method	Natural convection	Natural convection	Natural convection	Natural convection
Operating altitude, max.	2000 m	2000 m	2000 m	2000 m
Protection degree	IP20, Type 1	IP20, Type 1	IP20, Type 1	IP20, Type 1
Ambient temperature	10 – 40 °C	10 – 40 °C	10 – 40 °C	10 – 40 °C
Temperature: Storage	-10 – 60 °C	-10 – 60 °C	-10 – 60 °C	-10 – 60 °C
Emission sound pressure level (EN ISO 2151)	45 dB(A)	45 dB(A)	45 dB(A)	45 dB(A)
Weight	4.3 kg	4.3 kg	4.3 kg	4.3 kg

Tbl. 28: Technical data, MVP 030-3 DC

16.3 Substances in contact with media

Pump parts	Substances in contact with media
Housing cover	Aluminum alloy (AlMgSi)
Head cover	Aluminum alloy (AlMgSi)
Diaphragm clamping disk	Aluminum alloy (AlMgSi)
Diaphragm	PTFE
Valves	FKM
Inlet	Aluminum alloy (AlMgSi)
Outlet / silencer	Aluminum / silicone rubber
Threaded hose couplings	Anodized aluminum
Hose	PE

Tbl. 29: Materials that make contact with the process media

16.4 Dimensions

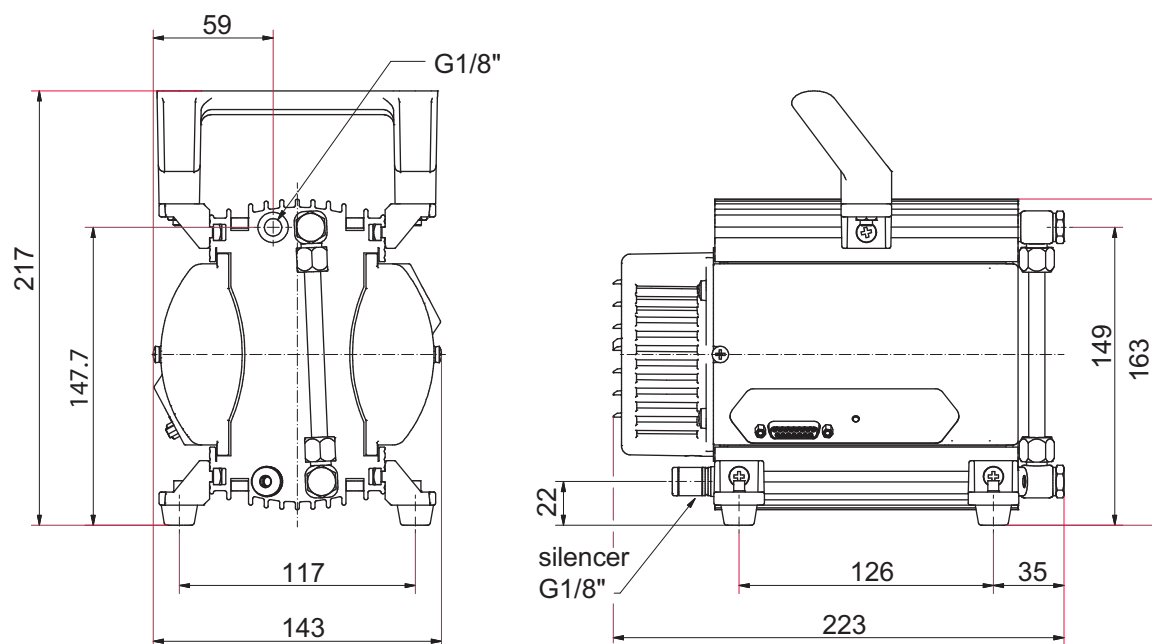


Fig. 15: Dimensions MVP 030-3 DC

Dimensions in mm

EC Declaration of Conformity

This declaration of conformity has been issued under the sole responsibility of the manufacturer.

Declaration for product(s) of the type:

Diaphragm pump

MVP 030-3 DC

We hereby declare that the listed product satisfies all relevant provisions of the following **European Directives**.

Machinery 2006/42/EC (Annex II, no. 1 A)

Electromagnetic compatibility 2014/30/EU

Restriction of the use of certain hazardous substances 2011/65/EU

Restriction of the use of certain hazardous substances, delegated directive 2015/863/EU

Harmonized standards and applied national standards and specifications:

DIN EN ISO 12100:2011

DIN EN IEC 63000:2019

DIN EN 1012-2:2011

DIN EN 61010-1:2020

DIN EN IEC 61326-1:2022

DIN EN ISO 80079-36:2016

DIN EN 1127-1:2019

The authorized representative for the compilation of technical documents is Dr. Adrian Wirth, Pfeiffer Vacuum GmbH, Berliner Strasse 43, 35614 Asslar, Germany.

Signature:



(Daniel Sälzer)
Managing Director

Pfeiffer Vacuum GmbH
Berliner Straße 43
35614 Asslar
Germany

Asslar, 2023-06-16



UK Declaration of Conformity

This declaration of conformity has been issued under the sole responsibility of the manufacturer.

Declaration for product(s) of the type:

Diaphragm pump

MVP 030-3 DC

We hereby declare that the listed product satisfies all relevant provisions of the following **British Directives**.

Supply of Machinery (Safety) Regulations 2008

Electromagnetic Compatibility Regulations 2016

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

Applied standards and specifications:

EN ISO 12100:2010

EN IEC 63000:2018

EN 1012-2:1996+A1:2009

EN 61010-1:2010 + A1:2019 + A1:2019/AC:2019

EN IEC 61326-1:2021

EN ISO 80079-36:2016

EN 1127-1:2019

The manufacturer's authorized representative in the United Kingdom and the authorized agent for compiling the technical documentation is Pfeiffer Vacuum Ltd, 16 Plover Close, Interchange Park, MK169PS Newport Pagnell.

Signature:



(Daniel Sälzer)
Managing Director

Pfeiffer Vacuum GmbH
Berliner Straße 43
35614 Asslar
Germany

Asslar, 2023-06-16

**UK
CA**

VACUUM SOLUTIONS FROM A SINGLE SOURCE

Pfeiffer Vacuum stands for innovative and custom vacuum solutions worldwide, technological perfection, competent advice and reliable service.

COMPLETE RANGE OF PRODUCTS

From a single component to complex systems:

We are the only supplier of vacuum technology that provides a complete product portfolio.

COMPETENCE IN THEORY AND PRACTICE

Benefit from our know-how and our portfolio of training opportunities!

We support you with your plant layout and provide first-class on-site service worldwide.

ed. H - Date 2511 - P/N: PU0076BEN



Are you looking for a
perfect vacuum solution?
Please contact us

Pfeiffer Vacuum GmbH
Headquarters • Germany
T +49 6441 802-0
info@pfeiffer-vacuum.de

www.pfeiffer-vacuum.com